



A12.1062

Laboratory Biological Microscope, Info LCD





A12.1062

Laboratory Biological Microscope, Info LCD

A12.1062 is the latest advanced laboratory microscope to be launched in 2021. It has a ergonomics and stable structure. The five hole encoded nosepiece, together with the base LCD screen, can provide such advanced functions as magnification display, brightness memory, and auto power off (ECO).

A12.1062's high-quality NIS45/60 series infinite flat field achromatic objective lens, super large 230x150mm working stage, 3W LED high brightness light source, and innovative 100x water lens enable high magnification observation to avoid the trouble of cleaning cedar oil.

Further more, A12.1062 has complete upgrade options (BF, DF, PH, PL, FL), making it competent for most laboratory microscopic observation work, it is an ideal choice for universities, laboratories and other fields.



Seidentopf Head

Inclined 30° Interpupillary Distance 47-78mm, Eyepiece Tube Dia.30mm

22mm Wide Field Of View

WF10x/22mm For A More Comprehensive Observation Content Sample Observation. The Eyepiece Adopts A Flat Field Distortion-free Design To Prevent The Edge Of The Field From Being Imaginary And Stray Light.



NIS60 Infinity Optical System

Nis Infinity Plan Objectives Can Provide High Contrast And Very Flat Image Up To FN 22. With WF0x/22mm Wide Field Eyepieces, The System Always Brings You Sharp, Excellent Resolution And High Signal To Noise Ratio Imaging.

A12.1062 Is Suitable For All Kinds Of Microscopic Using , Especially For Beginners And The Users With Long Time Micro-operation. The A12.1062 Series Of Microscopes Are Fully Optimized For The Needs Of Such Users. In Terms Of The Objectives, The Image Quality And Ease Of Use Are Achieved.

Plan Objective

By Using Infinity Plan Objectives, Flat Image With Higher Imaging Reduction Degree Over The Entire Field Of View Could Be Achieved.






100x Water-immersion Objective

Ordinary 100x Oil-immersion Objective Needs To Use Cedar Oil As The Observation Medium. After Use, It Needs To Be Cleaned With Ether Alcohol Or Xylene, Which Is Easy To Cause Air Pollution And Improper Cleaning. The Water - Immersion Objective Uses Water As A Medium To Solve The Above Problems Perfectly, Reducing The Damage To The Body And Environmental Pollution.

40x LWD Objective

The Working Distance Of 40x Objective Can Be Up To 1.5mm, Avoiding The Erosion From Residual Immersion Oil And Water When Converted From 100x To 40x Objective.

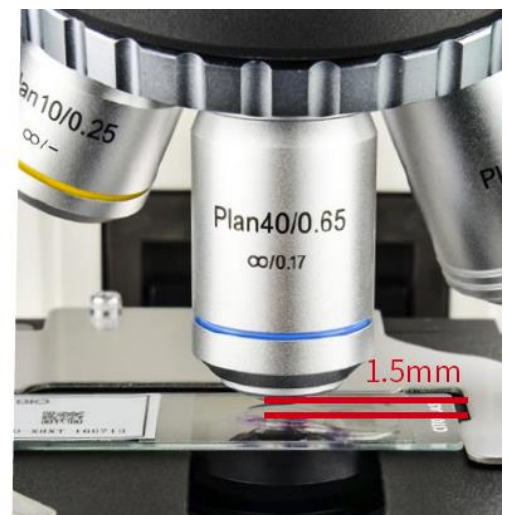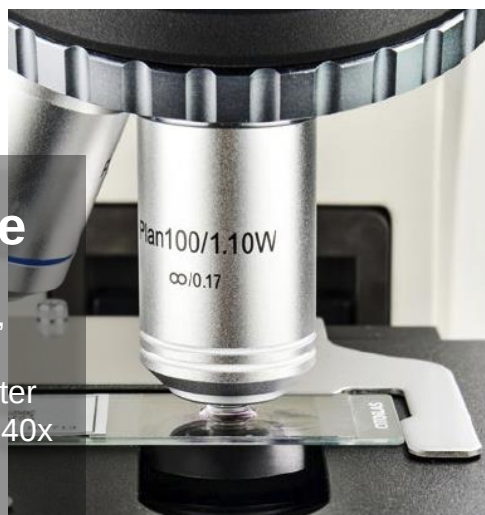

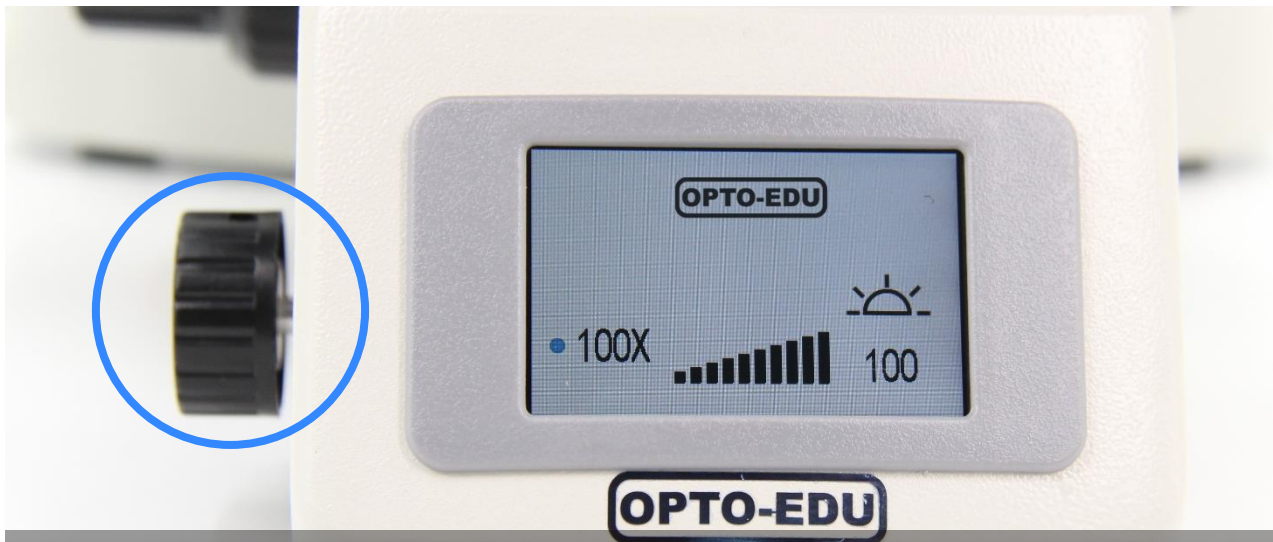


Use A Dimming Knob To Achieve Multiple Functions (Available In A12.1062 Series)


One Click: Enter Standby Status; Press + Up-spin: Switch To The Upper Light Source; Double Clicks: Light Lock Or Unlock; Rotation: Adjust Brightness; Press + Down-spin: Switch To The Under Light Source Press 3 Seconds: Set The Time Of Turning Off The Light After Leaving

The Display Of Microscope Use State (Available In A12.1062 Series)


The LCD Screen On The Front Of The Microscope Can Display The Using Status Of The Microscope, Including Magnification, Light Intensity, Sleepy Model, And So On.

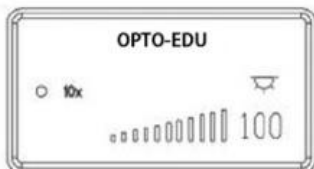


Start & Working Mode

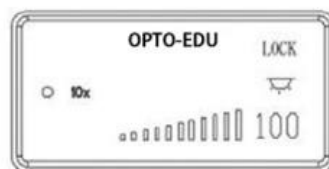

Lock Mode

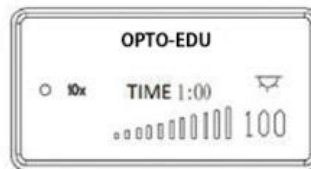

ECO Mode

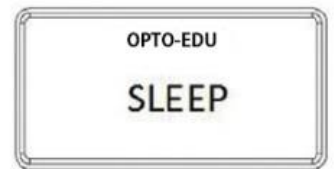

Sleep Mode




Coded Nosepiece (Available In A12.1062 Series)


It Can Memorize The Illumination Brightness When Using Each Objective. When Different Objectives Are Converted To Each Other, The Light Intensity Is Automatically Adjusted To Reduce Visual Fatigue And Improve Work Efficiency.




4x

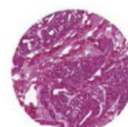


10x

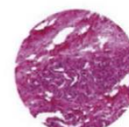

20x

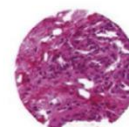

40x

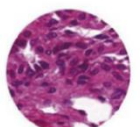

100x



Rackless Mechanical Working Stage

Double Layer Mechanical Rackless Stage, Size 230x150mm, Moving Range 78x54mm, Anti-Scratch Hard Coating Dural Platform



Abbe Condenser & Kohler Illumination

Inserted Abbe Condenser N.A. 1.25, With Slot For Phase Contrast Slide And Dark Field Slide, Work With 3W S-LED Kohler Illumination

A12.1061 Upgrade To A12.1062



A12.1061

A12.1062

A12.1061 Upgrade To A12.1062

- Objective NIS45 Upgrade To NIS60
- Condenser Upgrade To Coded Nosepiece
- Stand Upgrade To With LCD Screen
- 1W LED Upgrade To 3W LED.

Various Observation Methods

With The Deepening Of Research In Various Fields, A Single Observation Mode Can No Longer Meet The Daily Scientific Research Work. As A Continually Upgradable Microscope, A12.1062 Can Be Extended On Basic Models To Show A Variety Of Observation Capabilities.

Observation Methods	Bright Field	Dark Field	Phase Contrast	Fluorescent	Simple Polarizing
	•	•	•	•	•

Multifunctional Universal Condenser (For Bright Field, Dark Field And Phase Contrast)

A12.1062 Offers Universal Condensers For Bright Field, Dark Field And Phase Contrast. The Observation Methods Could Be Quickly Switched By Switching The Slider. The Phase Contrast And Bright Field Slider Is Universal For 4x- 100x Objectives Also, Simple And Fast To Use. The N.A. Value Index On The Aperture Diaphragm Of The Condenser Is Easily Set To Get Exact Size Of Diaphragm To Correspond With The Different Objectives.




LED EPI-Fluorescent Illumination

Can Upgrade To Fluorescent Microscope A16.1062. Led Epi - Fluorescent Illumination Is Safety And Convenient. You Don't Need Time To Warm Up Or Cool Down. You Don't Need To Align The Bulb, And The Long Lifetime Of Led Bulb Is Up To 5000 Hours. There Are Two Filters Position Available And Switching Is Fast And Easy.

Simple Polarizing Set

The installation of simple polarizing accessories is as shown in the figure. The polarizer is installed in the host lens barrel, and the polarizer is installed on the chassis light source. Slowly rotate the polarizer by hand to make it reach the orthogonal position of polarized light, and then the sample can be detected.

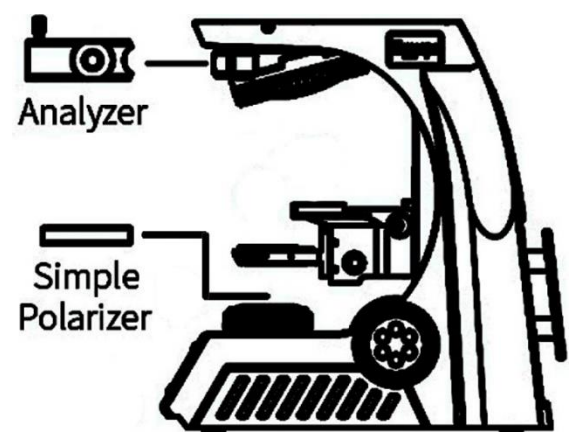




A12.1061, A12.1602 Laboratory Microscope A16.1062 LED Fluorescent Microscope		A12.1061		A12.1062		A16.1062		Cata. No.
		-B	-T	-B	-T	-B	-T	
Optical System	NIS Infinite Optical System	●	●	●	●	●	●	
Observation Method	Bright Field	●	●	●	●	●	●	
	Dark Field	○	○	○	○	○	○	
	Polarizing	○	○	○	○	○	○	
	Phase Contrast	-	-	○	○	○	○	
	Epi-Fluorescence	-	-	○	○	●	●	
Head	Seidentopf Binocular Head, Inclined At 30° , Interpupillary Distance 47-78mm, Eyepiece Tube Dia.30mm	●		●		●		A53.1061-B
	Seidentopf Trinocular Head, Inclined At 30° , Interpupillary Distance 47-78mm, Eyepiece Tube Dia.30mm, Light Path Fixed 50:50		●		●		●	A53.1061-T
	Seidentopf Trinocular Head, Inclined At 30° , Interpupillary Distance 47-78mm, Eyepiece Tube Dia.30mm, Light Path Switchable 0:100/100:0	○	○	○	○	○	○	A53.1061-TS
	Tilting Binocular Head, Inclined 5°~35° , Interpupillary Distance 47-78mm, Eyepiece Tube Dia.30mm	○	○	○	○	○	○	A53.1061-BS
	Digital Head, Built-in Digital Camera 5.0M, Support Android, IOS, Windows, Output Image By WIFI In Real Time	-	-	○	○	-	-	A53.1061-5.0M
	Upgrade to A31.1062-5.0M							
	Digital Head, Built-in Wireless Digital Camera 5.0M, Support Android, IOS, Windows, Output Image By WIFI In Real Time, App Software, By Scanning QR Code On Microscope, Installing App & Identifying Microscope, You Can View Micro Image On Mobile Phone & Tablet.	-	-	○	○	-	-	A53.1061-5.0W
Eyepiece	Upgrade to A31.1062-5.0W							
	EW10x/22mm, Diopter Adjustable, Dia.30mm	●	●	●	●	●	●	A51.1029-1022
	EW15x/16mm, Diopter Adjustable, Dia.30mm	○	○	○	○	○	○	A51.1029-1516
Nosepiece	EW20x/12mm, Diopter Adjustable, Dia.30mm	○	○	○	○	○	○	A51.1029-2012
	Quintuple Nosepiece, Dovetail Interface	●	●	-	-	-	-	
	Coded Quintuple Nosepiece, Dovetail Interface	-	-	●	●	●	●	

A12.1061, A12.1602 Laboratory Microscope A16.1062 LED Fluorescent Microscope		A12.1061		A12.1062		A16.1062		Cata. No.
		-B	-T	-B	-T	-B	-T	
NIS45 Infinity Plan Objective	4x/0.10, W.D.20.6mm	●	●	-	-	-	-	A52.1045-4
	10x/0.25, W.D.17.9mm	●	●	-	-	-	-	A52.1045-10
	20x/0.40, W.D.6.40mm	○	○	-	-	-	-	A52.1045-20
	40x/0.60, W.D.1.50mm	●	●	-	-	-	-	A52.1045-40
	60x/0.80, W.D.0.30mm	○	○	-	-	-	-	A52.1045-60
	100x/1.25(Oil), W.D.0.16mm	●	●	-	-	-	-	A52.1045-100
	100x/1.10(Water), W.D.0.16mm	○	○	-	-	-	-	A52.1045-100W
NIS60 Infinity Plan Objective	4x/0.10, W.D.30.0mm	-	-	●	●	●	●	A52.1060-4
	10x/0.25, W.D.10.2mm	-	-	●	●	●	●	A52.1060-10
	20x/0.40, W.D.4.8mm	-	-	○	○	○	○	A52.1060-20
	40x/0.65, W.D.1.5mm	-	-	●	●	●	●	A52.1060-40
	60x/0.80, W.D.0.30mm	-	-	○	○	○	○	A52.1060-60
	100x/1.25(Oil), W.D.0.3mm	-	-	●	●	●	●	A52.1060-100
	100x/1.10(Water), W.D.0.2mm	-	-	○	○	○	○	A52.1060-100W
NIS60 Infinity Plan Phase Contrast Objective	10x/0.25, W.D.10.2mm	-	-	○	○	○	○	A5C.1060-10
	20x/0.40, W.D.12.0mm	-	-	○	○	○	○	A5C.1060-20
	40x/0.65, W.D.0.7mm	-	-	○	○	○	○	A5C.1060-40
	100x/1.25(Oil), W.D.0.2mm	-	-	○	○	○	○	A5C.1060-100
Focusing	Coaxial Coarse & Fine Adjustment, Tension Adjustable, Fine Division 0.002mm, Coarse Focusing Range 28mm	●	●	●	●	●	●	
Working Stage	Double Layer Mechanical Rackless Stage, Size 230x150mm, Moving Range 78x54mm	●	●	-	-	-	-	
	Double Layer Mechanical Rackless Stage, Size 230x150mm, Moving Range 78x54mm, Anti-Scratch Hard Coating Dural Platform	-	-	●	●	●	●	
Condenser	Inserted Abbe Condenser N.A. 1.25, With Slot For Phase Contrast Slide And Dark Field Slide.	●	●	●	●	●	●	
Illumination	1W S-LED Critical Illumination	●	●	-	-	-	-	
	3W S-LED Kohler Illumination	-	-	●	●	●	●	
	Green Filter	●	●	●	●	●	●	
Dark Field	Dark Field Slide	○	○	○	○	○	○	A5D.1062
Polarizing	Simple Polarizing Set	○	○	○	○	○	○	A5P.1008
Phase Contrast	Phase Slider For 10x-40x	○	○	○	○	○	○	A5C.1062-S1040
	Phase Slider For 100x	○	○	○	○	○	○	A5C.1062-S100
Epi-Fluorescence Attachment	3W LED Epi-Fluorescence Attachment, Fly-Eye Lens Illumination, With 2 Filter Position + Bright Field Position, Switch Quickly, With Attachable UV Shield, B, G Filters	○	○	○	○	●	●	A5F.1062
	Filter U	○	○	○	○	○	○	A5F.1062-U
	Filter V	○	○	○	○	○	○	A5F.1062-V
Photo Adapter	1.0x C-Mount	○	○	○	○	○	○	A55.1062-1.0
	0.5x C-Mount	○	○	○	○	○	○	A55.1062-0.5
Power Supply	AC 100-240V,50/60Hz	●	●	●	●	●	●	
	USB Charging Port, Support To Use Exeternal Rechargeable Battery Power Bank Like Mobile Phone Out Door Freely	○	○	○	○	○	○	A56.1060-U
LCD Screen	LCD Screen On Front Of Body, Display Using State Of Microscope, Including Magnification, Light Intensity, Standby Status, Set Power Off Timer 5 Mins to 8 Hours, And So On.	-	-	●	●	●	●	
	Light Color Temperature Adjustable 3000K-7000K	-	-	○	○	○	○	A56.1060-C
Dimensions	220(W)x290(D)x472(H) mm	●	●	●	●	●	●	

Note: "●" In Table Is Standard Outfits, "○" Is Optional Accessories "-" Is Unavailable

A12.1061 Has The Multifunctional Digital Head, The User Does Not Have To Be Confined In Front Of The Microscope. Instead, It Can Be Used For Mobile Microscope Teaching And Outdoor Field Observation Through Mobile Terminals And External Mobile Power. The Objective, Eyepiece, And Observation Tube Are Effectively Anti-mold Treated, So You Can Ensure A Consistently Clear Image And Extend The Life Of The Microscope, Even When Working In Hot And Humid Environments.

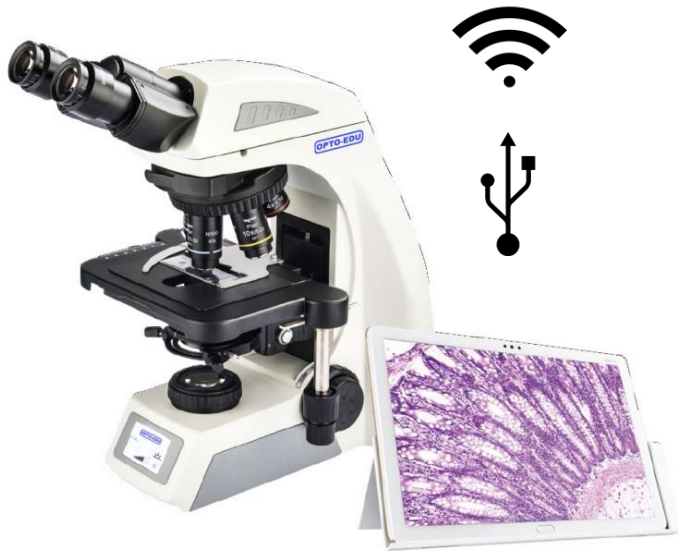

A31.1062 Digital Microscope, WIFI & USB

Built-in Camera 5.0M, USB 2.0 or WIFI Output, Supporting Android, IOS, Windows Operating System, Wired And Wifi Modes The Image Under The Microscope Can Be Output To The External Device In Real Time, And There Is No Data Line Connection, And The Operator Can Move More Freely.

Professional Microscopic Imaging Software

Microscopic Imaging Observation, Analysis And Processing Can Be Performed On External Devices, Including Photographing, Measurement, Image Adjustment, Storage, Synthesis, Etc.



Mobile Devices Perform Image Browsing And Processing By Scanning Code

A31.1062-5.0W WIFI Digital Microscope, By Scanning The Qr Code On The Microscope, Installing The App And Identifying The Microscope, You Can View The Microscopic Image On Your Phone And Tablet.



External Rechargeable Battery



A USB Charging Port Is Reserved On The Body, Which Can Be Used As A Microscope Power Source. This Microscope Can Also Be Used Outside And During Power Outages To Get Rid Of The Microscope's Dependence On The Power Outlet.






The Microscope Is Compact And Can Be Placed In An Ordinary Classroom Closet. It Has A Special Carrying Handle, And Is Also Lightweight And Stable. The Microscope Back Plate Is Designed With A Hub Device To Effectively Store The Long Power Cord, Improve The Cleanliness Of The Laboratory, And Reduce The Tripping Accident Caused By The Long Power Cord During The Carrying Process. The Wooden Storage Box Is An Optional Accessory That Is Very Convenient For Storage And Carrying.

Ergonomic Design





In Daily Scientific Research Teaching And Pathological Diagnosis, Working In Front Of The Microscope For A Long Time Has Become The Norm, And The Consequent Use Fatigue Often Leads To Physical Discomfort, Thereby Reducing Work Efficiency And Effectiveness. This A12.1061 Microscope Uses An Ergonomic Design, High Eye-point, Low-hand Focus Mechanism, Low-hand Stage And Other Ergonomic Designs To Ensure The User Can Perform Microscope Operation In The Most Comfortable Situation. . The Focus Knob, Illumination Control Knob And Stage Handle Are All In Close Proximity. The User Can Put Both Hands On The Table While Working, And Can Operate A12.1061 With Minimal Movement.


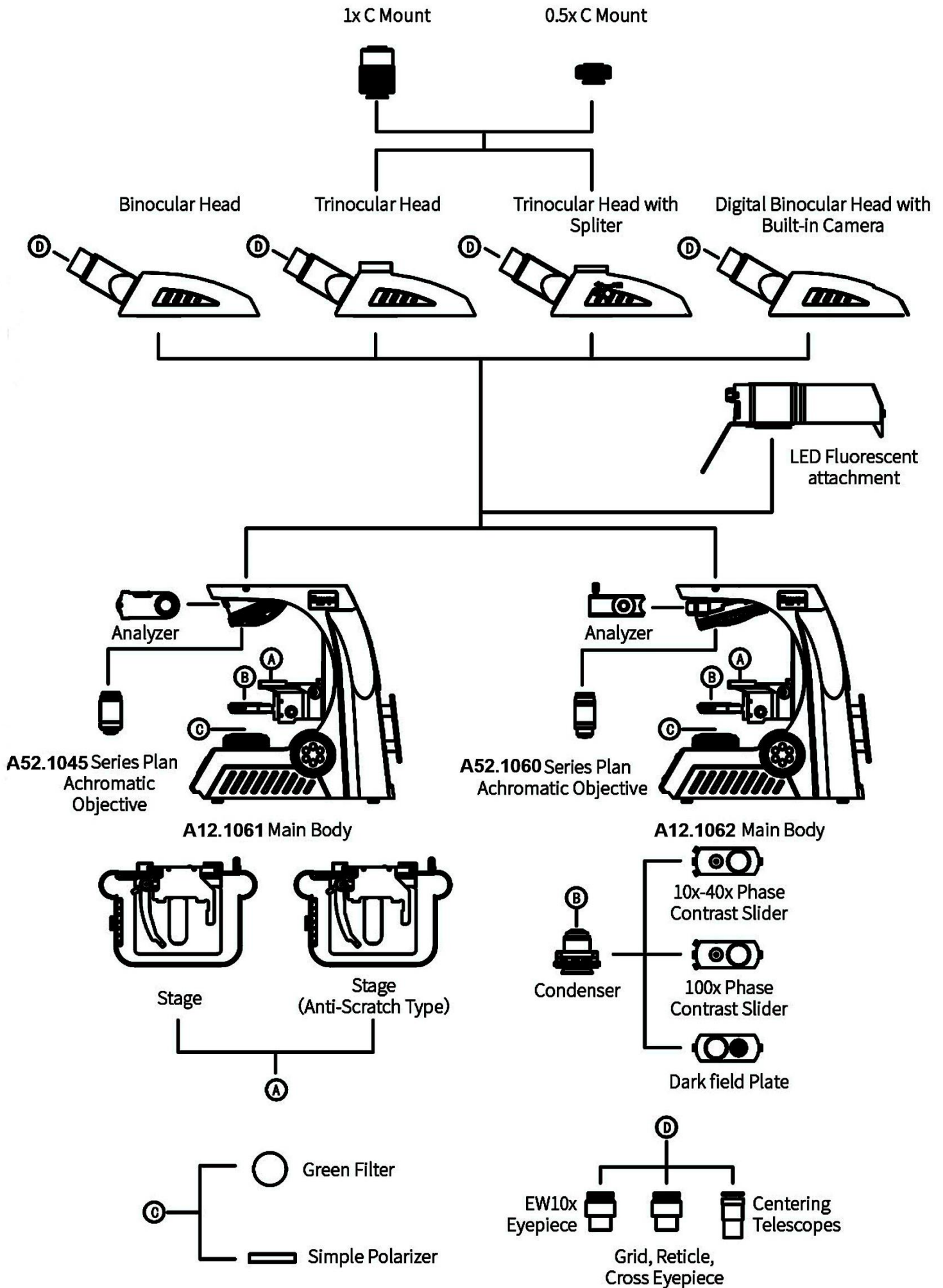


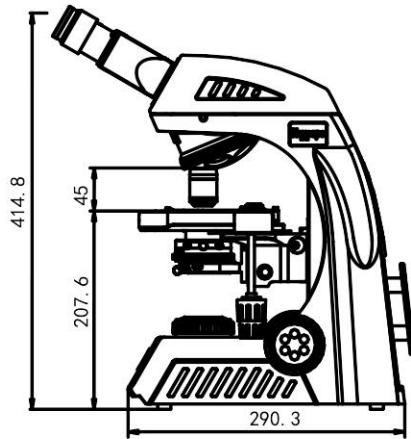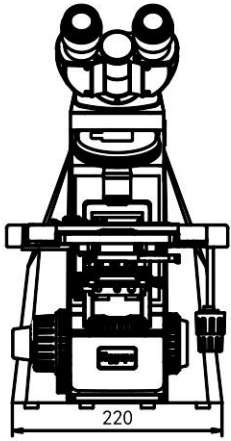


A12.1061

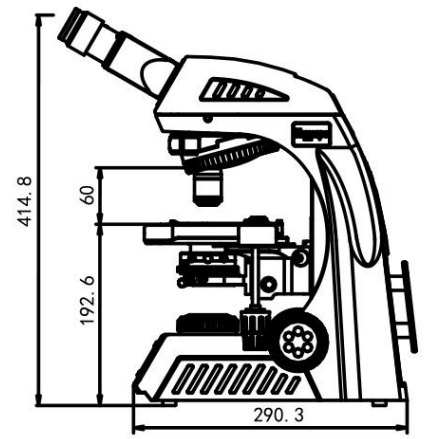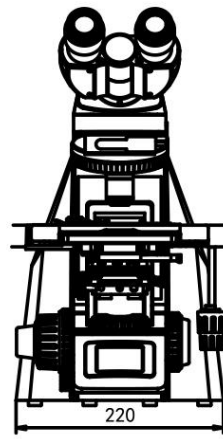

A12.1062

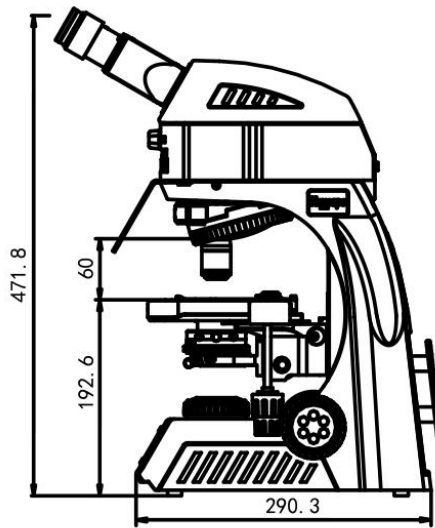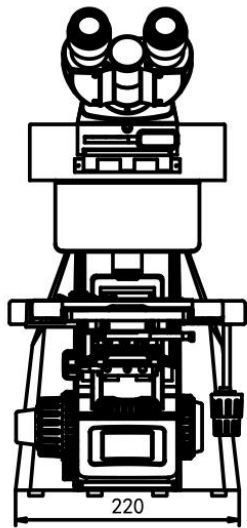


A16.1062

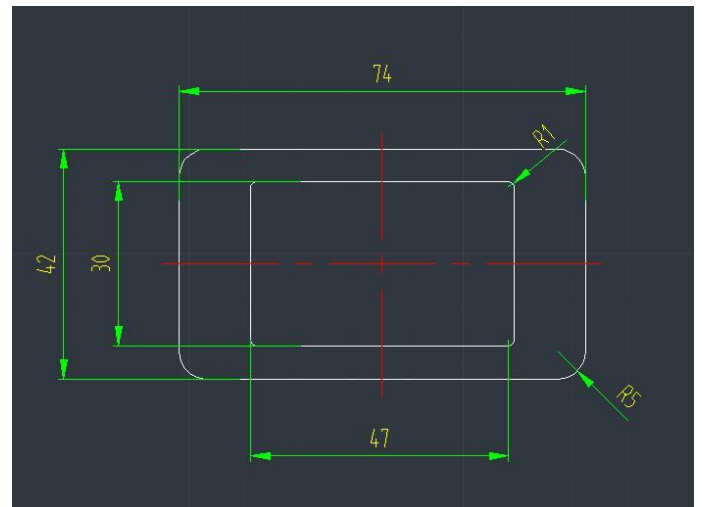

Info LCD

A12.1062 Packing



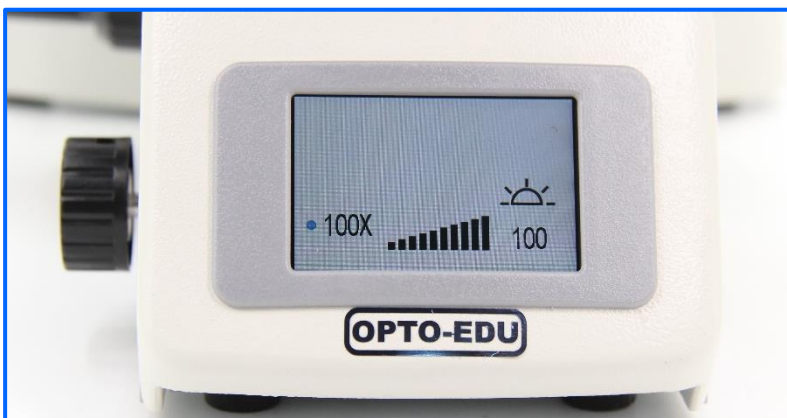

MICROSCOPE IS OUR FOCUS

The Key To Micro World

Opto-Edu is one of the most professional supplier for microscope & educational instruments from China. We have been focusing in this field for more than 25 years, has a professional sales team who can assist our customer to find the BEST SELECTED microscope with competitive price and 3 YEARS Warranty.

25+ Year professional experience we know Microscopes the best!

150+ Microscope & accessory manufacturers supply all models from China

200+ Hot sale microscopes & Newest Models Updated Every Month

750+ Customer from all over the world & keep rising every day

1500+ Microscope products create your one-stop purchase platform

3000+ Educational Instruments For School, College And University Teaching

The Most Professional Microscope Manufacturer in China!



MICRO
WORLD

The Key To OPTO-EDU



Opto-Edu (Beijing) Co., Ltd.

F-1501 Wanda Plaza, No.18 Shijingshan Road, Beijing 100043, China

Official Main Website: www.optoedu.com

www.cnoec.com.cn, www.optoedumicroscope.com, www.microscopemadeinchina.com

Skype: xincnoec Wechat, Mobile, WhatsApp: +86 13911110627

Tel: +86 10 88696085 Email: sale@optoedu.com

